

## AUTOMATED ELECTROPHORESIS

# Experion™ Pro260 Analysis Kit

## Quick Guide

### 1 Temperature Equilibration

Remove the Pro260 stain (**blue cap**), Pro260 sample buffer (**yellow cap**), Pro260 ladder (**red cap**), and 2 tubes of Pro260 gel (**green cap**) from storage. Allow these reagents to warm to room temperature (~15 min). Briefly vortex and then centrifuge the tubes before use to collect the solutions at the bottom of the tubes.

### 2 Prepare the Gel and Gel-Stain Solution

- To prepare the gel-stain solution, add 20 µl Pro260 stain (**blue cap**) to a tube of Pro260 gel (520 µl; **green cap**); vortex briefly
- Pipet the gel-stain solution and the contents of the other tube of Pro260 gel (520 µl; **green cap**) into separate spin filters; centrifuge at 10,000 x g for 5 min

Note: Unused filtered gel-stain solution and filtered gel may be stored for up to 1 month at 4°C in the dark. Unused gel should be refiltered after 1 month.

### 3 Prepare the Sample Buffer

#### Reducing Conditions

- Add 1 µl 2-mercaptoethanol to 30 µl sample buffer (**yellow cap**) for each chip to be run; prepare fresh solution daily

#### Nonreducing Conditions


- Add 1 µl deionized water (0.2 µm filtered) to 30 µl sample buffer (**yellow cap**) for each chip to be run. In addition, prepare a tube of sample buffer under reducing conditions as instructed above; the Pro260 ladder always requires a reducing sample buffer

### 4 Prepare the Samples and Pro260 Ladder

- Add 2 µl sample buffer to 4 µl each sample; vortex
- Add 2 µl sample buffer to 4 µl Pro260 ladder (**red cap**); vortex
- Heat the samples and ladder at 95–100°C for 3–5 min; centrifuge the samples briefly
- Add 84 µl deionized water (0.2 µm filtered) to samples and ladder; vortex

### 5 Prime the Chip

- Remove a Pro260 chip from its packaging and place it on the chip platform in the Experion priming station
- Add 12 µl gel-stain solution to the gel priming well
- Close the lid of the priming station, set the pressure to **B**, and set the time to **3**
- Press the **Start** button
- When priming is complete, remove the primed chip from the priming station
- Flip the chip over, and visually inspect the microchannels for trapped air bubbles or incomplete priming

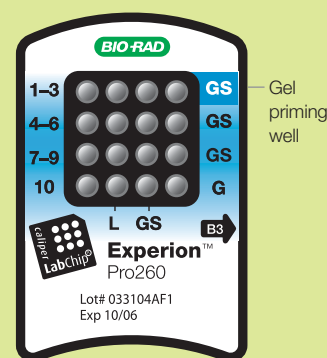



# Experience Meets Innovation

- 6 Load the Samples and Pro260 Ladder Into the Chip**
- Remove and discard any remaining gel-stain solution from the gel priming well
  - Pipet 12  $\mu$ l filtered gel-stain solution into all 4 wells labeled **GS**
  - Pipet 12  $\mu$ l filtered gel into the well labeled **G**
  - Pipet 6  $\mu$ l diluted protein ladder into the well labeled **L**
  - Pipet 6  $\mu$ l diluted sample into each sample well. Do not leave any sample well empty. If necessary, pipet a sample or ladder replicate into any empty sample wells
  - Run the chip in the Experion electrophoresis station within 5 min of loading

- 7 Run the Pro260 Analysis**
- Turn the Experion electrophoresis station on and then launch the Experion software
  - Place the chip on the chip platform of the electrophoresis station and close the lid
  - Select **New Run**, then **Experion Protein 260** Assay
  - Click the **Start** button on the screen to begin the run
  - When the run is complete, remove and discard the used chip

- 8 Clean the Electrodes**
- Place a cleaning chip filled with 800  $\mu$ l deionized water (0.2  $\mu$ m filtered) into the electrophoresis station
  - Close the lid for 60 sec to clean the electrodes, and then open it for 60 sec to allow the electrodes to dry completely
  - Remove the cleaning chip

## Ordering Information

Catalog #	Description
700-7000	Experion System, 100–120/220–240 V, for protein analysis, includes electrophoresis station, priming station, software, USB2 cable, test chip, instructions (analysis kits sold separately)
700-7101	Experion Pro260 Analysis Kit for 10 Chips
700-7102	Experion Pro260 Analysis Kit for 25 Chips
700-7151	Experion Pro260 Chips, 10
700-7152	Experion Pro260 Reagents and Supplies, for 10 chips
700-7251	Experion Cleaning Chips, 10
700-7254	Experion Spin Filters, 10
163-2091	ReadyPrep™ Proteomics Grade Water, 500 ml
161-0710	2-Mercaptoethanol, 25 ml
700-7256	Experion Pro260 Ladder



LabChip and the LabChip logo are trademarks of Caliper Life Sciences, Inc. Bio-Rad Laboratories, Inc. is licensed by Caliper Life Sciences, Inc. to sell products using the LabChip technology for research use only.

These products are licensed under US Patent Nos. 5,863,753, 5,658,751, 5,436,134, and 5,582,947, and pending patent applications, and related foreign patents, for internal research and development use only in detecting, quantitating, and sizing macromolecules, in combination with microfluidics, where internal research and development use expressly excludes the use of this product for providing medical, diagnostic, or any other testing, analysis, or screening services, or providing clinical information or clinical analysis, in any event in return for compensation by an unrelated party.